



BMA BIOMEDICALS

BMA BIOMEDICALS
Rheinstrasse 28-32
CH-4302 Augst (Switzerland)
Phone:++41 61 811 6222
Fax: ++41 61 811 6006
info@bma.ch
www.bma.ch

Royalmount Aqueous Mounting Medium

Royalmount is an aqueous mounting medium for cover-slipping tissue sections stained with alcohol soluble chromogens like fast red or AEC (3-amino-9-ethylcarbazole).

Product Number:	T-3402/250ml
Quantity:	250ml
	Supplied as liquid solution ready to use
Stability:	Original vial: 1 year at room temperature
Applications:	Before using Royalmount remove the red cap of the bottle and cut off the tip of the nozzle with a sharp blade or a scissors. Make sure that the opening is not too wide. Do not shake the bottle to avoid air bubbles.

Typical Protocol:

- Remove excess liquid from the slide. Tissue must still be humid.
- Apply 1-3 drops of Royalmount to the tissue section.
- Place slide horizontally on a heating plate at 60°C for at least 30 minutes.
- Remove slide from the heating plate and allow it to reach room temperature.
- To remove Royalmount, soak slide overnight in water.
- If required, post-mounting a coverslip on the hardened Royalmount is possible. Apply a few drops of a commercial mounting medium for coverslips on the slide and mount the coverslip.

Royalmount is also available in 60ml size (product T-3402/60ml)

Please see www.bma.ch for protocols and general information.

For in vitro research only.

T-3402/250ml

Royalmount

10.11.2011